LBX ELFV – 750 – 001

Please read the User Manual carefully before use and follow all operating and safety instructions

user manual

english, español

User Manual

LBX ELFV

Preface

Thank you for purchasing our product. Users should read this manual carefully, follow the instructions and procedures, and beware of all the preventive measures when using this instrument.

Service

If help is needed, you can always contact your dealer or Labbox via www.labbox.com

Please, provide the customer service representative with the following information:

- Serial number
- Description of the problem
- Your contact information

Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 12 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts that have been damaged due to improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

SAFETY PRECAUTIONS

WHEN USED CORRECTLY, THESE UNITS POSE NO HEALTH RISK.

HOWEVER, THESE UNITS CAN DELIVER DANGEROUS LEVELS OF ELECTRICITY AND ARE TO BE OPERATED ONLY BY QUALIFIED PERSONNEL FOLLOWING THE GUIDELINES LAID OUT IN THIS INSTRUCTION MANUAL.

ANYONE INTENDING TO USE THIS EQUIPMENT SHOULD READ THE COMPLETE MANUAL THOROUGHLY.

THE UNIT MUST NEVER BE USED WITHOUT THE SAFETY LID CORRECTLY IN POSITION.

THE UNIT SHOULD NOT BE USED IF THERE IS ANY SIGN OF DAMAGE TO THE EXTERNAL TANK OR LID.

Cleaning and maintenance

HORIZONTAL UNITS

- Use warm water, but not above 60°C, as it can damage the unit and components
- Use a mild detergent or water with a mild concentration of soap.
- Do not leave the units in detergent for more than 30 min.
- Rinse the tank thoroughly with warm water or distilled water to prevent buildup of salts.
- Be careful not to damage the enclosed electrode.
- Vigorous cleaning is neither necessary nor advised.
- Let the components air dry before use.
- Compatible detergents include dishwashing liquid, Hexane and Aliphatic hydrocarbons.
- **Never let the following cleaning agents come into contact with the unit:**
 - Acetone
 - Phenol
 - Chloroform
 - Carbon tetrachloride
 - Methanol
 - Ethanol
 - Isopropyl alcohol

RNase DECONTAMINATION

This can be done by following the instructions below:

- Clean the units with mild detergent as described above
- Wash with 3% hydrogen peroxide (H₂O₂) for 10 min.
- Rinse with 0.1% DEPC (diethyl pyrocarbonate) treated with distilled water.

CAUTION: DEPC is a suspected carcinogen. Always take the necessary precautions when using it.

Operation instructions


SETTING UP THE VERTICAL GEL TANKS

Instructions for fitting the Electrode Cables:


1. Check the position of the lid on the unit. This shows the polarity and the correct orientation of the cables: black is negative and red is positive.
2. Remove the lid from the unit. Note that if the lid is not removed, fitting the cables may cause the gold plug to loosen and damage the electrode.
3. Screw the cables into the tapped holes as tightly as possible so that there is no gap between the lid and the leading edge of the cable fitting.
4. Refit the lid.

VERTICAL GEL CASTING

1. Clean the glass plates – rinse them with distilled water and let them air dry.
2. Overlap a spacer glass with a notched one.
3. Put the two plates in the core. Keep their bottoms perfectly aligned and in contact with the table.
4. Insert a cuneal plate closely outside the two plates.
5. Do the same for the other side.
Note: the spacer plate must be placed in the outer side and the notched one in the inner side.
6. Pull out the handle of the gel casting base and rotate it downwards.
7. Put the core inside the gel casting base.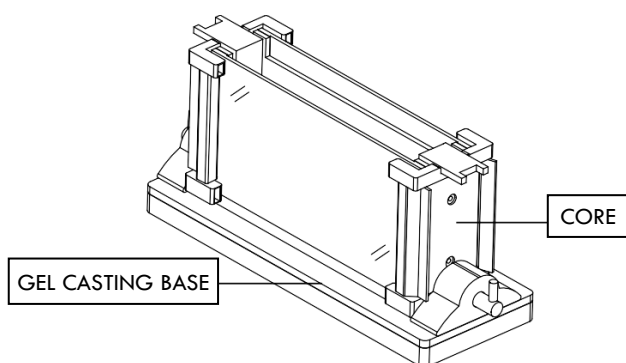

8. Push the handle into the whole and rotate upwards.

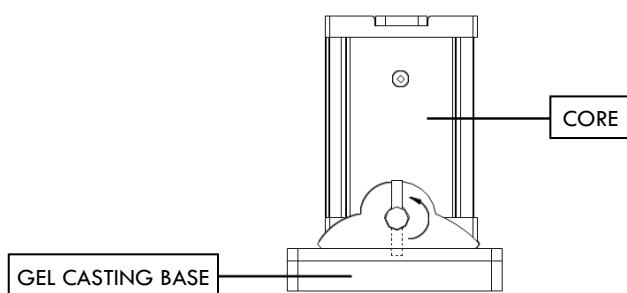GEL POURING

1. Pour the gel into the blocks. Do it carefully to avoid generating bubbles.
2. Place the comb(s) carefully. Make sure there are no air bubbles under the ends of the comb teeth.
3. Let the gel settle and wait for it to set completely.
4. Pull the comb(s) out carefully and move the tray with the gel into the main tank.

RUNNING THE GEL

1. Put the core inside the bottom tank.
2. Fill both tanks with 1 x buffer solution.
3. Load the samples into the wells using pipettes. Be careful not to damage the wells or induce bubbles.
4. Carefully cover the tank with the lid and connect it to a voltage supply.
5. Run the electrophoresis. Note that gels are typically run under 150-200V. Higher voltages allow for a faster electrophoresis but with poorer quality results.

Warranty

This product has a one-year warranty that covers any defects in the material and workmanship the replacement of any part (except the platinum wire) should any defect come up during the warranty period. However, the warranty is voided under the following circumstances:

- Defects caused by improper operation.
- Repair or modification done by anyone other than the supplier or an authorized agent.
- Use with cables or connectors not specified by the supplier.
- Deliberate or accidental misuse.
- Damage caused by disaster.

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