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Please read the User Manual carefully before use, and follow all operating and safety instructions!

user manual

english / español / français

User Manual

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Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

Service

In order to guarantee this equipment Works safely and efficiently, it must have a regular maintenance. In case of any faults, do not try to repair it yourself. If help is needed, you can always contact your dealer or Labbox via www.labbox.com

Please, provide the customer service representative with the following information:

- Serial number
- Description of the problem
- Your contact information

Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

Warning: do not use the rotor if it has been damaged due to corrosion and/or fatigue.

1. Safety Precautions

The precautions described in the User Manual are carefully developed with intent to cover all the possible risks. Please read them carefully before operating the centrifuge.

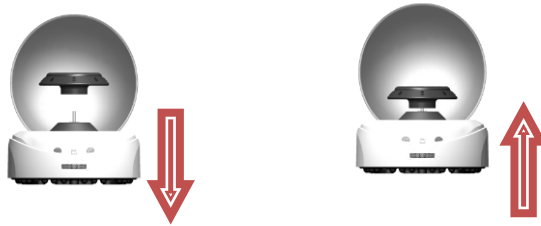
- Ensure that only trained staff works with the instrument.
- Never use the centrifuge in any way other than that specified in these instructions manual.
- Never place your hands in the rotor area unless the rotor has completely stopped.

If something abnormal occurs during the operation of this centrifuge, please stop it immediately and contact your supplier/manufacturer.

2. Loading the Rotor

Line up the rotor with the motor shaft. Press the rotor down to secure it with the shaft.

- Please use your both hands to load the rotor.

3. Operation

- Open the lid and press the ON/OFF switch located at the back of the unit.
- Symmetrically load the centrifuge tubes on the rotor.
- Close the centrifuge's lid to begin running.
- Switch off, the rotor will gradually stop.
- Open the lid and do not remove the samples until the rotor completely stops moving.

4. Specifications

Max. Speed [rpm]	7000
Max. Relative centrifugal force [x g]	2680
Rotor Capacity	1) 8×0.5/1.5/2.0ml Tubes 2) 16 x 0.2ml PCR tubes 2×0.2ml PCR 8-tube Strip
Run Time	Continuous
Driving Motor	DC motor
Power	AC 100V~240V, 50Hz/60Hz
Noise Level	≤45 dB
Dimensions [D×W×H mm]	160 x 170 x 122
Weight [kg]	0.5

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